

Vent Kit Assembly Details

Code 663-100 to 663-108

Australia

1800 358 101
sales.au@westlab.com

Canada

1 877 822 1455
sales.ca@westlab.com

New Zealand


09 972 7682
sales.nz@westlab.com

United States

646 568 5391
sales.us@westlab.com

Installation

1. Screw nipple & reducer into vent opening at base of cabinet.
2. Slide weight support onto reducer ensuring weight is transferred to floor, rather than cabinet.
3. Screw elbow onto reducer.
4. Place vent pipe through roof & ceiling penetration & screw into elbow.
5. Cut off vent pipe from 300mm to 1000mm min height above roof.
6. Rivet adaptor to vent pipe.
7. Rivet whirlybird to adaptor.
8. Seal proof penetration.

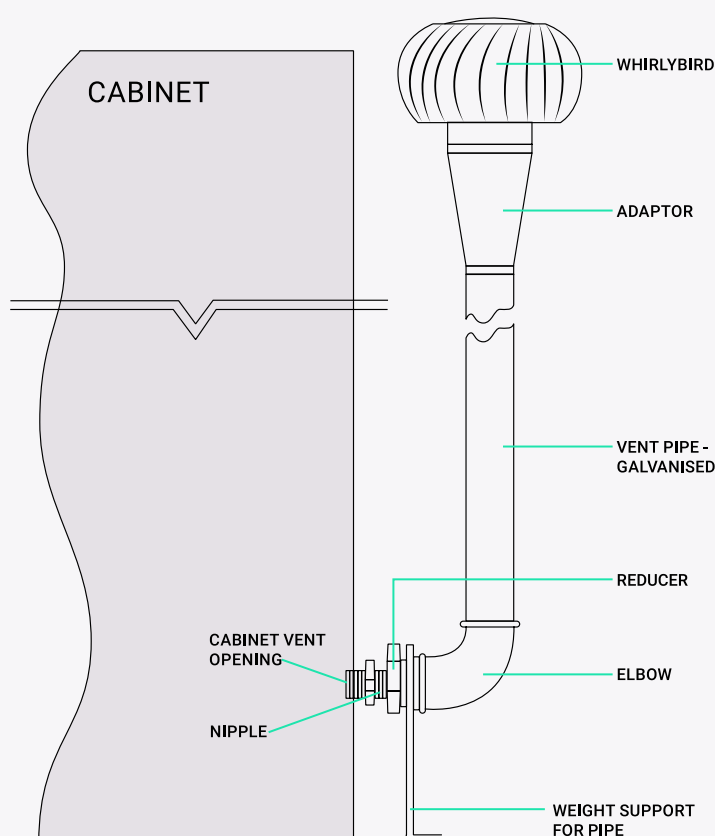

VENT ASSEMBLY KIT DETAILS

**CODE 663-100
to 663-108**

ORDERING WITH WESTLAB IS EASY

Online

www.westlab.com.au

Order online with our easy Express Order form for fast, secure ordering.

Phone

1800 385 101

Speak to one of our friendly, efficient staff from

8.00am - 5.00pm (Mon/Fri)

Email

sales@westlab.com.au

24 hours, 7 days a week convenience.